
Human Cytokine Magnetic 10-Plex Panel Technical Data Sheet

Performance Characteristics of the Human Cytokine Magnetic 10-Plex Panel (Cat. No. LHC0001M)

Typical Standard Curves as generated for Lot 1 (Note: Data presented here were collected as a 10-plex, but are depicted in two 5-plex groupings for clarity.)

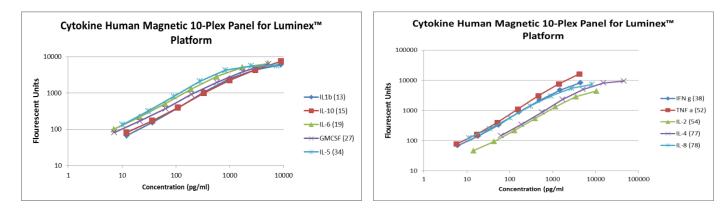

| | IL-1β | IL-10 | IL-6 | GM CSF | IL-5 | IFN γ | TNF α | IL-2 | IL-4 | IL-8 |
|---------------------------|----------|----------|----------|---------|-----------|-----------|-----------|---------|----------|----------|
| Magnetic Bead Region | 13 | 15 | 19 | 27 | 34 | 38 | 52 | 54 | 77 | 78 |
| Sensitivity | < 1 | < 0.5 | < 0.5 | < 0.5 | < 0.5 | < 0.5 | < 0.5 | < 1 | < 1 | < 1 |
| Recovery (Serum) | ٧ | v | v | v | v | ٧ | ٧ | v | v | v |
| Recovery (Heparin Plasma) | ٧ | ٧ | ٧ | ٧ | v | NR | ٧ | v | v | ٧ |
| Recovery (EDTA Plasma) | ٧ | v | v | v | v | ٧ | v | v | v | v |
| Recovery (Citrate Plasma) | ٧ | v | ٧ | v | v | ٧ | ٧ | v | v | ٧ |
| NIBSC calibration | 1ng=25IU | 1ng=13IU | 1ng=30IU | 1ng=6IU | 1ng=0.2IU | 1ng=0.7IU | 1ng=0.2IU | 1ng=7IU | 1ng=65IU | 1ng=40IU |
| NIBSC Preparation Number | 86/680 | 93/722 | 89/548 | 88/646 | 90/586 | 82/587 | 88/786 | 86/564 | 88/656 | 89/520 |
| Inter assay variation | 8.8% | 9.8% | 7% | 9.1% | 7.5% | 9% | 8.3% | 9.5% | 8.7% | 9.8% |

Notes: 70-130% recovery (v), 50-69% recovery (-), 131-150% recovery (+), <50% or >150% recovery (NR-not recommended)

Limited product warranty

Life Technologies Corporation and/or its affiliate(s) warrant their products as set forth in the Life Technologies' General Terms and Conditions of Sale found on Life Technologies' website at **www.thermofisher.com/us/en/home/global/terms-and-conditions.html**. If you have any questions, please contact Life Technologies at **www.thermofisher.com/support**.



Manufacturer's address: Bender MedSystems GmbH | Campus Vienna Biocenter 2 | 1030 Vienna, Austria

The information in this guide is subject to change without notice.

DISCLAIMER: TO THE EXTENT ALLOWED BY LAW, LIFE TECHNOLOGIES AND/OR ITS AFFILIATE(S) WILL NOT BE LIABLE FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE, OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING YOUR USE OF IT.

Important Licensing Information: These products may be covered by one or more Limited Use Label Licenses. By use of these products, you accept the terms and conditions of all applicable Limited Use Label Licenses.

©2022 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified.

For Research Use Only. Not for use in diagnostic procedures. MAN0017197 Rev. 2.0 (01) Effective Date: 31 January 2022 Manufacturing site: Bender MedSystems GmbH | Campus Vienna Biocenter 2 | 1030 Vienna, Austria **Thermo Fisher** S C I E N T I F I C www.thermofisher.com